

Dissolution of Sodium Hydroxide Pellets using a LevMixer™

Mixing system: Newmix®- Levtech® LevMixer™

Mixing bag: 100L L-Mix bag

Mixing type: Powder-liquid

The Newmix-Levtech LevMixer is a compact and non-invasive single-use mixing system. The heart of this system is a mixing bag incorporating a bottom-mounted levitating impeller designed for powder-liquid and liquid-liquid mixing applications. The impeller is frictionless and generates no particles. All product-contacting surfaces are 100% disposable.

Introduction

Powder-liquid mixing is a common requirement in biopharma processing. In order to maximize mixing speed for powder-liquid applications, various impeller sizes and locations are available.

In this experiment, a LevMixer was used to prepare 70L of a buffer solution typically encountered in bioprocess facilities.



Hoegaarden, Belgium - Europe

Phone: +32 (0) 16.76.61.59

Lexington, KY - USA

Phone: 859.263.1135

Minneapolis, MN - USA

Phone: 952.942.0855

www.atmi-lifesciences.com

info@atmi-lifesciences.com

© 2008 ATMI, Inc. All Rights Reserved.

ATMI, the ATMI logo, LevMixer, Levtech, the LevTech logo and Newmix are trademarks or registered trademarks of Advanced Technology Materials, Inc in the United States, other countries or both.

Application note XA208E 0807rev1

Experimental

A 100L LevMixer mixing bag, fitted with a large (6.35") centrally-mounted impeller, was filled with 70L of water, and mixing speed was set to 214rpm. Once a vortex had formed, 2.5kg of sodium hydroxide (NaOH) pellets was added to the mixing bag, resulting in an approximately 35g/L solution.

Results

Dissolution was observed visually. After 5 minutes, all of the NaOH pellets were seen to have dissolved. The temperature of the water increased from initial 30°C to 38°C during this experiment.

Conclusions

The Newmix-Levtech LevMixer system is well suited to preparation of buffer solutions such as sodium hydroxide. A powder-liquid mixing bag, which includes a centrally-mounted, large, levitated impeller, is a good choice for such applications. Mixing times in the sub-300 second range are typical at the tested volume and concentration.

In view of the temperature increase observed, ATMI recommends that the impeller be started prior to adding the NaOH; this will reduce the chance of local hotspots that could damage the bag material. The use of a jacketed tank and chiller can be considered to if rapid cooling is desirable.